

Analytical and Testing Instruments for the HPI Market

Shimadzu's Total Solutions for the Hydrocarbon Processing Industry












World Map of Shimadzu Sales, Service, Manufacturing, and R&D Facilities



Shimadzu's Total Solutions for the Hydrocarbon Processing Industry (Petrochemical, Chemical)

The Hydrocarbon Processing Industry (HPI), encompassing petroleum refining, gas processing, petrochemical and chemical, is a foundational field for all industry. In the HPI market, many kinds of analytical and measuring instruments are used for quality control purposes and process management.

With a wide range of products, from GC and Elemental Analyzers to an On-line Water Quality Analyzer, Shimadzu offers total support for laboratories in the HPI market. From Upstream to Downstream, our rugged, high-quality products are perfectly suited to meet the demands of your laboratory and contribute to your business success.

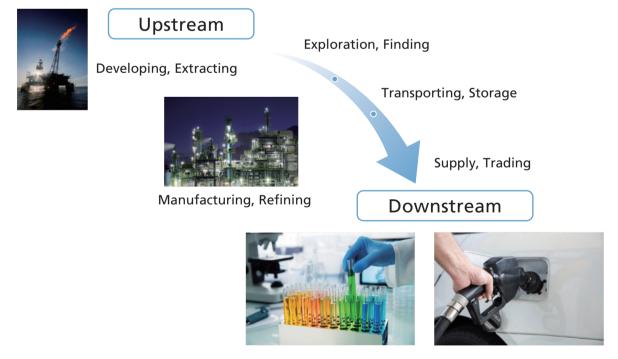



For more information about Shimadzu's solutions for the HPI market, visit our website at: https://www.shimadzu.com/an/industry/petrochemicalchemical/index.html

Shimadzu Customized GC Analyzers

Various Applications in the HPI Market

- Support third-party detectors, including a pulsed flame photometric detector (PFPD) and a pulsed discharge helium ionization detector (PDHID).
- Support third-party software for detailed hydrocarbon analysis (DHA) and Chrom Merge.
- Support liquefied petroleum gas (LPG) sampling devices and vaporizers.
- Complete application manual and methods



Nexis GC-2030

Product Lineup

	Natural Gas Analyzers	Refinery Gas Analyzer (RGA)	Process and Custom Gas Analyzers	TOGA/DGA Analyzer	Petrochemicals analyzer
Applicable Methods	ASTM D1945/D1946 GPA 2261, GPA 2177 GPA 2286, GPA 2186 ISO 6975	UOP 539, ASTM D7833, D 2163	ASTM D1946 ASTM E1746	ASTM D3612C	ASTM D6228
Analysis	C1-C6+ Hydrocarbons, Non Condensable gases O2, N2, CO, CO2 and H2S Optional Hydrogen Extended NGA C1-C14+ NGL, Dual Oven NGA High Speed NGA	C1-C6+ Hydrocarbons saturates and unsaturates H2S, Non Condensable gases O2, N2, CO, CO2 and H2 LPG Sampling Vaporizer	O2, N2, CO, CO2 Impurities in Chlorine	TOGA with Head Space H2, O2, N2, CO, CO2, C1, C2's, C3's	Trace Sulfur
Features	Analysis time 10 min.	Analysis time less than 6 min.	Special Hastelloy Material for Chlorine Analyzer	High Sensitivity Extension up to C4	PFPD for Sulfur

Transformer Oil Gas Analyzer (TOGA)



Simulated Distillation Analyzer




Trace Sulfur Analysis GC System

Sulfur Chemiluminescence Detection Gas Chromatograph System

Nexis SCD-2030


- Best-in-Class Sensitivity
- Unmatched Stability performance
- Excellent Maintainability
- Automation Functions

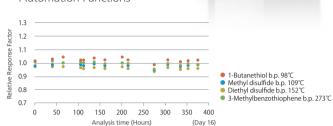

Excellent Long-Term Stability

VUV PIONA+ Analysis GC System

- Comprehensive gasoline measurements determining bulk PIONA and specific oxygenate and BTEX compound content.
- Provides a lot of information in a fast, robust measurement.
- Relatively simple instrumentation, setup, and analysis procedures.

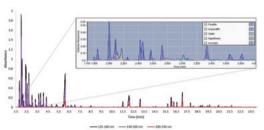

VUV Chromatogram of Gasoline Sample


* The VUV detector VGA-100 is a product of VUV ANALYTICS. INC.



Simulated Distillation	DHA Analyzers	Oxygenates Analyzers	Sulfur Analyzers	Aromatics Analyzers
ASTM D7096, D7213, D3710, D2887, D6417, D6352, D7500, D5307, D7169, IP 480, ISO3924	ASTM D5134, D6729, D6730, D6733, D8071	ASTM D4815,D7423 UOP 960	ASTM D6228 ASTM D5504 ASTM D5623 ASTM D7011	ASTM D3606, 5580
Boiling Point Distribution C1-C110 (As per method)	C1-C14 Hydrocarbons PIONA Report	C1-C4 Alcohols	Trace Sulfur Analysis Individual components	Aromatics
Fast Simdist Correlations with D86 and D1160, Integrated Shimadzu Software	Tuned D6730 column Fast DHA Front End PIONA Room Temperature Database VUV PIONA+	Use of N2 Carrier for D4815	Use of FPD, PFPD or SCD Detectors	Combination model for D3606, D5580. D4815

Trace Moisture Analysis GC System

- Easily analyzes both liquid and gas samples
- Uses a variety of Watercol[™] GC columns depending on the matrices
- Uses a simple, high-sensitivity, and very stable low-cost Barrier Ionization Discharge (BID) Detector

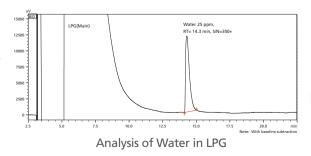



Aromatics Analysis HPLC



Features

Switching Valve Configuration

Super Performance Expandability


Precise Sample Injection

Applicable Methods

ASTM D 6591, D6379, IP 391, IP 436, IP 548, IP590 Custom Methods

Analysis Subjects

Aromatics in fuels Poly-aromatics in diesel FAME in jet fuel

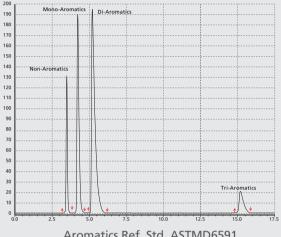

Aromatics Ref. Std. ASTMD6591


Applicable Methods

ASTM D5837

Custom Methods

Analysis Subjects

Components in transformer oil

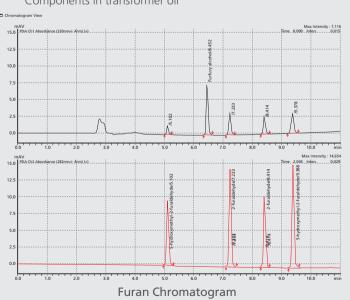

Furan Analysis HPLC



LC/MS/MS



LCMS-8060

Features




Scan Speed 30000u / sec



- Applicable Methods
 - Custom Methods
- Analysis Subjects

Additives in polymers

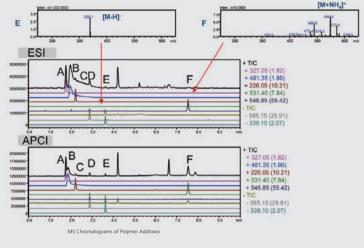

Analysis of Additives in Plastic

Applicable Methods EPA 8270B, TO -14, IP-585

Custom Methods

Analysis Subjects

Volatile organic compounds Arsine phosphine in ethylene FAME in jet fuel



HS-20 Headspace Analysis System



TD-30 Thermal Desorption System

VOCs Analysis GC-MS



GCMS-QP2020 NX

Features



Inductively Coupled Plasma Mass Spectrometer (ICP-MS)



Features






High Stability, High Sensitivity and Low Interference

Applicable Methods

Customized

Analysis Subjects

Trace elements in naphtha Determination of toxic arsenic species using LC-ICP-MS

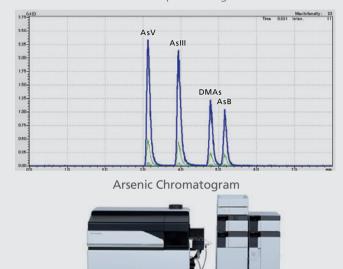


Applicable Methods

Elemental Analysis ASTM D4951, D5184 ASTM D4951,D5184 UOP 389, UOP 303 UOP 714, APHA 3120, IFP 9507

Analysis Subjects

All elements Lubricating oils Petroleum naphtha, gasoline Marine fuel, fuel oil, heavy oils

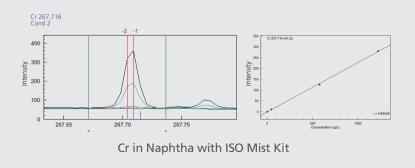

ICP Atomic Emission Spectrometer (ICPE)





Atomic Absorption Spectrophotometer (AAS)



Features

Unique 3D Double Beam Optics




Dual Background Correction Functions

Applicable Methods

ASTM D6595, D3237, D3567, D3605, D3635, D3831, D3919, D4628, IFP9312, UOP946, D5863

Analysis Subjects

Additives in polymers All elements Lubricating oil Petroleum oils Lead in gasoline

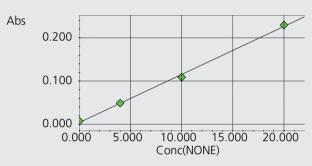


Abs=0.011100Conc+0.0043463 r=0.9989

Arsenic in Vacuum Gas Oil (VGO)

- Applicable Methods ASTM D5839, ASTM D4294, C114 ISO 8754
- Analysis Subjects

Sulfur in oil
All elements

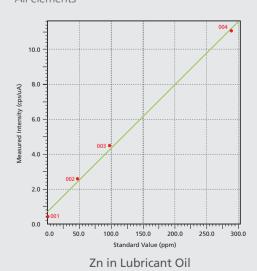

Energy Dispersive X-ray Fluorescence Spectrometer (EDX)



EDX-7000/8000/8100

Features







Fourier Transform Infrared Spectrophotometer (FTIR)



Features






- Applicable Methods
 D7371, EN 14078, D7575, D7678

 Custom Methods
- Analysis Subjects

Fame in biodiesel Total petroleum hydrocarbons Additives in polymers Oil in water

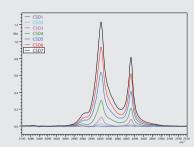



Enlarged View of Peaks Around 2920 cm⁻¹

Calibration Curve and Standard Sample Concentrations


Analysis of Oil in Water


Applicable Methods ASTM D1840, D2008

Custom Methods

Analysis Subjects

Petroleum products Turbine fuels Mineral oils

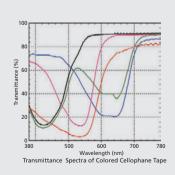

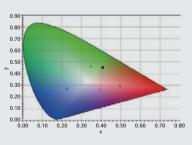

Chromaticity Diagram of XY Color System

Analysis of Colored Cellophane Tape

UV-VIS Spectrophotometer (UV-Vis)



Features





High Performance to Meet Diverse Needs

Thermal Analyzers



DSC-60 Plus

Features



Diverse Measurements by Simple Operations




Applicable Methods

ASTM D3417, D3418, D4419, D4591, D5483, ASTM D3850, D6370, D6375, **Custom Methods**

Analysis Subjects

Thermal properties of polymers DSC, TGA, TG/DTA, TMA

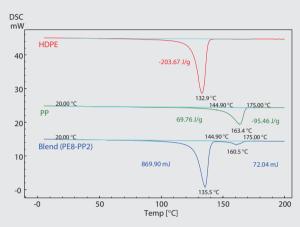


DSC Curves of HDPE, PP, and Blend Sample

Applicable Methods


ASTM D5412, JPI-5S-71-2010 Custom Methods

Analysis Subjects

Measuring oil in water Coumarin in diesel

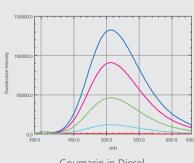

Coumarin in Diesel

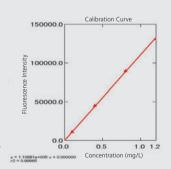

Spectrofluorophotometer



Features





Efficiency

Total Organic Carbon (TOC) Analyzer




Features



- Applicable Methods
 - ASTM D7573, EN 1484, EPA 415.1 APHA 5310B
- Analysis Subjects

TOC in various types of water, including high salt content water

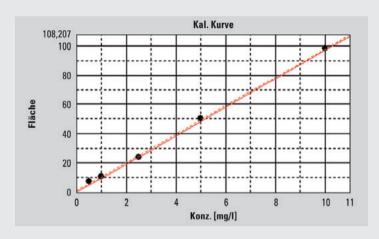


TOC in Brine Solution

Applicable Methods

Custom Methods

Analysis Subjects

TOC in various types of water

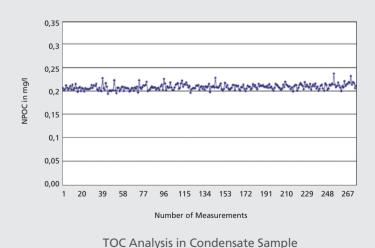

Mean Value 0.208mg/L, SD 0.006mg/L, DL 0.05mg/L

On-line TOC Analyzer



Features



Low Operation Costs



Testing Machines




AG-Xplus Series

Features




Eight Fold High Resolution

- Applicable Methods ISO 6892, ISO 527, ASTM D638
- **Analysis Subjects** Tensile properties of metals Tensile properties of plastics

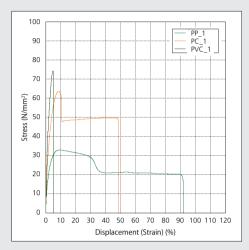

Tensile Test of PP, PC, PVC

Applicable Methods

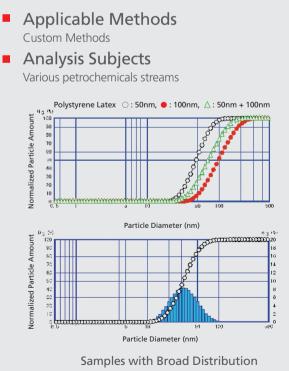

Particle Size Analyzer



SALD-2300

Features








Shimadzu Analytical Network Data System

LabSolutions CS


Controlling other vendor's GC and HPLC using LabSolutions DB/CS

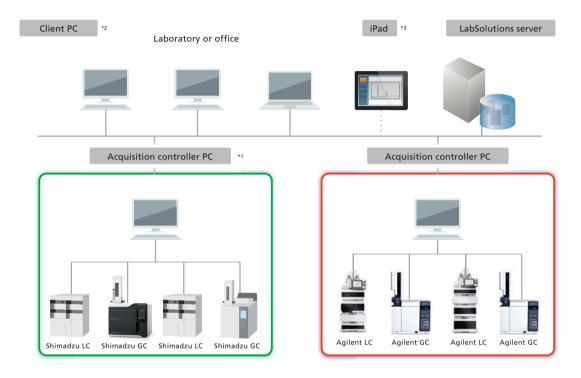

- *1 The acquisition controller PC controls the analytical instruments.

 Analysis directions and re-analysis of data can be performed using a client PC.
- *2 It is not necessary to install LabSolutions software on the client PC for terminal service.
- ¹³ When using an iPad, the installation of Citrix's XenApp is required. iPad is a registered trademark of Apple Inc.

Simple User Interface

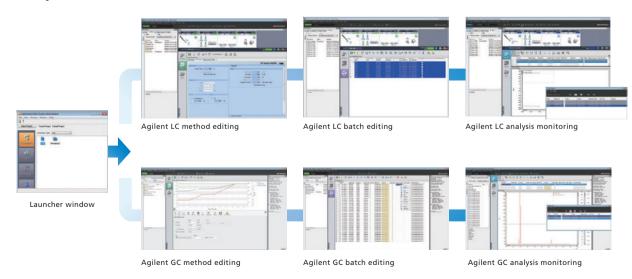



Sulfur Chemiluminescence Detection Gas Chromatograph System

Nexis SCD-2030

The Next Industry Standard SCD

The Nexis SCD-2030 is a next-generation sulfur chemiluminescence detection system. It has been developed to fulfill the unmet needs of laboratories the world over. The dramatically enhanced sensitivity and reliability, the excellent maintainability, and the automation functions, a first for the industry, will improve laboratory productivity.

- ► The New Standard in Reliability
- Dramatically Improved Productivity
- ▶ Best-in-Class Sensitivity



Shimadzu Corporation www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

Company names, products/service names and logos used in this publication are trademarks and trade names of Shimadzu Corporation, its subsidiaries or its affiliates, whether or not they are used with trademark symbol "TM" or "®".

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The contents of this publication are provided to you "as is" without warranty of any kind, and are subject to change without notice. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication.