

Application Data Sheet

No. 173

System Gas Chromatograph

Detailed Hydrocarbon Analysis of Naphtha Nexis GC-2030PONA2

A representative sample of the Naphtha is introduced into a gas chromatogram equipped with a methyl silicone bonded phase fused silica capillary column. Helium carrier gas transports the vaporized sample through the column in which the components are separated. Components are sensed by a FID as they elute from the column. Each eluting peak is identified and by comparing its retention index to a table of retention indices and by visual matching with a standard chromatogram. The mass concentration of each component is determined by area normalization with response factors. Peaks eluting after n-nonane are summed and reported as C₁₀⁺.

Analyzer Information

System Configuration:

One SPL Injector / one capillary column / one FID

Sample Information:

Methane
Ethane
Propane
Isobutane
n-Butane
Neopentane
Isopentane
n-Pentane
2,2-Dimethylbutane
Cyclopentane
2,3-Dimethylbutane
2-Methylpentane
3-Methylpentane
n-Hexane
2,2-Dimethylpentane
Methylcyclopentane
2,4-Dimethylpentane
2,2,3-Trimethylbutane
Benzene
3,3-dimethylpentane
Cyclohexane
2-Methylhexane
2,3-Dimethylpentane
1,1-Dimethylcyclopentane
3-Methylhexane

Concentration Range:

No.	Name of Compound	Concentration Range	
		Low Conc.	High Conc.
1	Methane	100 ppm	-
2	Ethane	100 ppm	-
3	Propane	100 ppm	-
4	Isobutane	100 ppm	-
5	n-Butane	100 ppm	-
6	Neopentane	100 ppm	-
7	Isopentane	100 ppm	-
8	n-Pentane	100 ppm	-
9	2,2-Dimethylbutane	100 ppm	-
10	Cyclopentane	100 ppm	-
11	2,3-Dimethylbutane	100 ppm	-
12	2-Methylpentane	100 ppm	-
13	3-Methylpentane	100 ppm	-
14	n-Hexane	100 ppm	-
15	2,2-Dimethylpentane	100 ppm	-
16	Methylcyclopentane	100 ppm	-
17	2,4-Dimethylpentane	100 ppm	-
18	2,2,3-Trimethylbutane	100 ppm	-
19	Benzene	100 ppm	-
20	3,3-dimethylpentane	100 ppm	-
21	Cyclohexane	100 ppm	-
22	2-Methylhexane	100 ppm	-
23	2,3-Dimethylpentane	100 ppm	-
24	1,1-Dimethylcyclopentane	100 ppm	-
25	3-Methylhexane	100 ppm	-

Sample Information:

cis-1,3-Dimethylcyclopentane
trans-1,3-Dimethylcyclopentane
3-Ethylpentane
trans-1,2-Dimethylcyclopentane
2,2,4-Trimethylpentane
n-Heptane
Methylcyclohexane + cis-1,2-Dimethylcyclopentane
1,1,3-Trimethylcyclopentane + 2,2-Dimethylhexane
Ethylcyclopentane
2,5-Dimethylhexane + 2,2,3-Trimethylpentane
2,4-Dimethylhexane
1,trans-2,cis-4-Trimethylcyclopentane
3,3-Dimethylhexane
1,trans-2,cis-3-Trimethylcyclopentane
2,3,4-Trimethylpentane
Toluene + 2,3,3-Trimethylpentane
1,1,2-Trimethylcyclopentane
2,3-Dimethylhexane
2-Methyl-3-ethylpentane
2-Methylheptane
4-Methylheptane + 3-Methyl-3-ethylpentane
3,4-Dimethylhexane
1,cis-2,trans-4-Trimethylcyclopentane + 1,cis-2,cis-4-Trimethylcyclopentane
cis-1,3-Dimethylcyclohexane
3-Methylheptane + 1,cis-2,trans-3-Trimethylcyclopentane
3-Ethylhexane + trans-1,4-Dimethylcyclohexane
1,1-Dimethylcyclohexane
2,2,5-Trimethylhexane + trans-1,3-Ethylmethylcyclopentane
cis-1,3-Ethylmethylcyclopentane

No.	Name of Compound	Concentration Range	
		Low Conc.	High Conc.
26	cis-1,3-Dimethylcyclopentane	100 ppm	-
27	trans-1,3-Dimethylcyclopentane	100 ppm	-
28	3-Ethylpentane	100 ppm	-
29	trans-1,2-Dimethylcyclopentane	100 ppm	-
30	2,2,4-Trimethylpentane	100 ppm	-
31	n-Heptane	100 ppm	-
32	Methylcyclohexane + cis-1,2-Dimethylcyclopentane	100 ppm	-
33	1,1,3-Trimethylcyclopentane + 2,2-Dimethylhexane	100 ppm	-
34	Ethylcyclopentane	100 ppm	-
35	2,5-Dimethylhexane + 2,2,3-Trimethylpentane	100 ppm	-
36	2,4-Dimethylhexane	100 ppm	-
37	1,trans-2,cis-4-Trimethylcyclopentane	100 ppm	-
38	3,3-Dimethylhexane	100 ppm	-
39	1,trans-2,cis-3-Trimethylcyclopentane	100 ppm	-
40	2,3,4-Trimethylpentane	100 ppm	-
41	Toluene + 2,3,3-Trimethylpentane	100 ppm	-
42	1,1,2-Trimethylcyclopentane	100 ppm	-
43	2,3-Dimethylhexane	100 ppm	-
44	2-Methyl-3-ethylpentane	100 ppm	-
45	2-Methylheptane	100 ppm	-
46	4-Methylheptane + 3-Methyl-3-ethylpentane	100 ppm	-
47	3,4-Dimethylhexane	100 ppm	-
48	1,cis-2,trans-4-Trimethylcyclopentane + 1,cis-2,cis-4-Trimethylcyclopentane	100 ppm	-
49	cis-1,3-Dimethylcyclohexane	100 ppm	-
50	3-Methylheptane + 1,cis-2,trans-3-Trimethylcyclopentane	100 ppm	-
51	3-Ethylhexane + trans-1,4-Dimethylcyclohexane	100 ppm	-
52	1,1-Dimethylcyclohexane	100 ppm	-
53	2,2,5-Trimethylhexane + trans-1,3-Ethylmethylcyclopentane	100 ppm	-
54	cis-1,3-Ethylmethylcyclopentane	100 ppm	-

Sample Information:

trans-1,2-Ethylmethylcyclopentane
2,2,4-Trimethylhexane +
1,1-Ethylmethylcyclopentane
trans-1,2-Dimethylcyclohexane
1,cis-2,cis-3-Trimethylcyclopentane
trans-1,3-Dimethylcyclohexane + cis-1,4-Dimethylcyclohexane
n-Octane
Isopropylcyclopentane + 2,4,4-Trimethylhexane
cis-1,2-Ethylmethylcyclopentane + 2,3,5-Trimethylhexane
2,2-Dimethylheptane
cis-1,2-Dimethylcyclohexane
2,2,3-Trimethylhexane
2,4-Dimethylheptane
4,4-Dimethylheptane
Ethylcyclohexane + n-Propylcyclopentane
2-Methyl-4-ethylhexane
2,6-Dimethylheptane
1,1,3-Trimethylcyclohexane
2,5-Dimethylheptane + 9P
3,5-Dimethylheptane + 3,3-Dimethylheptane
Ethylbenzene
Unidentified Naphthene + 2,3,4-Trimethylhexane
m-Xylene
p-Xylene
2,3-Dimethylheptane
3,4-Dimethylheptane + N
3,4-Dimethylheptane
4-Ethylheptane + N
4-Methyloctane
2-Methyloctane
3-Ethylheptane + N

Methods met:
ASTM-D5134

No.	Name of Compound	Concentration Range	
		Low Conc.	High Conc.
55	trans-1,2-Ethylmethylcyclopentane	100 ppm	-
56	2,2,4-Trimethylhexane + 1,1-Ethylmethylcyclopentane	100 ppm	-
57	trans-1,2-Dimethylcyclohexane	100 ppm	-
58	1,cis-2,cis-3-Trimethylcyclopentane	100 ppm	-
59	trans-1,3-Dimethylcyclohexane + cis-1,4-Dimethylcyclohexane	100 ppm	-
60	n-Octane	100 ppm	-
61	Isopropylcyclopentane + 2,4,4-Trimethylhexane	100 ppm	-
62	Ethylmethylcyclopentane + 2,3,5-Trimethylhexane	100 ppm	-
63	2,2-Dimethylheptane	100 ppm	-
64	cis-1,2-Dimethylcyclohexane	100 ppm	-
65	2,2,3-Trimethylhexane	100 ppm	-
66	2,4-Dimethylheptane	100 ppm	-
67	4,4-Dimethylheptane	100 ppm	-
68	Ethylcyclohexane + n-Propylcyclopentane	100 ppm	-
69	2-Methyl-4-ethylhexane	100 ppm	-
70	2,6-Dimethylheptane	100 ppm	-
71	1,1,3-Trimethylcyclohexane	100 ppm	-
72	2,5-Dimethylheptane + 9P	100 ppm	-
73	3,5-Dimethylheptane + 3,3-Dimethylheptane	100 ppm	-
74	Ethylbenzene	100 ppm	-
75	Unidentified Naphthene + 2,3,4-Trimethylhexane	100 ppm	-
76	m-Xylene	100 ppm	-
77	p-Xylene	100 ppm	-
78	2,3-Dimethylheptane	100 ppm	-
79	3,4-Dimethylheptane + N	100 ppm	-
80	3,4-Dimethylheptane	100 ppm	-
81	4-Ethylheptane + N	100 ppm	-
82	4-Methyloctane	100 ppm	-
83	2-Methyloctane	100 ppm	-
84	3-Ethylheptane + N	100 ppm	-

Detection limits may vary depending on the sample.
Please contact us for more consultation.

System Features

- Single FID channel
- Good repeatability
- PIONA Report by Dragon DHA Software*

Typical Chromatograms

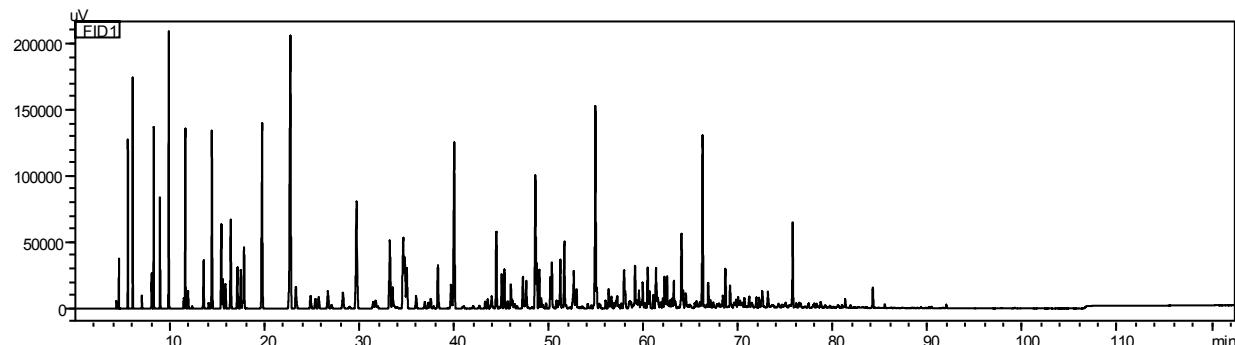


Fig. 1 Chromatogram of FID

SUMMARY REPORT			
Group Type	Total(Mass%)	Total(Vol%)	Total(Mol%)
Paraffins:	16.09	17.27	19.12
I-Paraffins:	19.99	21.17	22.47
Olefins:	0.10	0.10	0.14
Naphthenes:	25.26	24.50	30.14
Aromatics:	8.90	8.38	9.92
Total C10+:	28.92	27.84	17.48
Total Unknowns:	0.75	0.73	0.73

Oxygenates:

Total:	0.00(Mass%)	0.00(Vol%)
Total Oxygen Content:	0.00(Mass%)	
Multisubstituted Aromatics:	4.22(Mass%)	3.73(Vol%)

Average Molecular Weight: 119.92
Relative Density: 0.73
Reid Vapor Pressure @ 100F: 2.10psi - 14.46kPa
Calculated Octane Number: 46.5
Motor Octane Number (Jenkins Calculation): 45.4

	IBP	T10	T50	T90	FBP
BP by Mass (Deg F)	82.11	174.54	258.22	488.66	488.66
BP by Vol (Deg F)	82.11	174.54	258.22	488.66	488.66

Percent Carbon: 85.54 Percent Hydrogen: 14.46
Bromine Number (Calc): 0.15

Fig. 2 Example of PIONA Report *

* Dragon DHA software is registered trademark of Envantage Inc.

First Edition: November,2017



Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.
The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

© Shimadzu Corporation, 2017